


KAPS Adhesive Sealer Cover Removal Procedure

1. Isolate KAPS sealer from power sources, Electricity and Pneumatic.
2. Remove transparent Foil Cover from Right side and retain.
3. Remove 6 cover retaining fixings, 3 fixings left side- 2 fixings right side and 1 fixing right side under foil cover as picture.



KAPS Adhesive Sealer Cover Removal Procedure

4. Lift cover vertically ½ way, to expose electrical connection of touch screen, care should be taken to avoid damage to the emergency stop button.
5. Unplug the touch screen Ribbon cable and protective earth lead (green) from PCB and chassis connections (pull off connections) see picture.

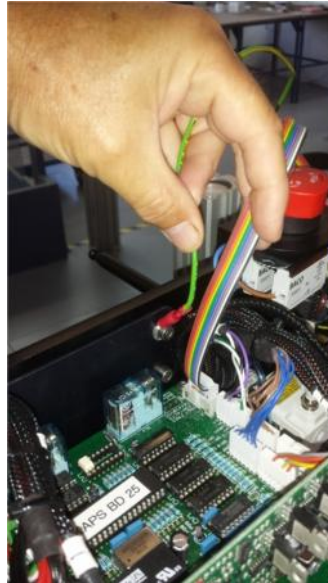

6. After removal of touch screen and protective earth connections, continue to lift cover vertically to fully remove cover from KAPs chassis.

