

Scorpion
Cover removal procedure

1. Disconnect all Power and Pneumatic services to Scorpion Plate stacker.
2. Remove Plate Hotels from Scorpion.
 - a) Slide Hotel Locking Bar to release one side of Plate Hotel.
 - b) Tilt Plate Hotel to remove.



3. Remove both hotel locking bars from Chassis (two screws each side)



4. Remove the six cover retaining screws.(three each side)



Scorpion
Cover removal procedure

5. Move gripper arm away from home position, to enable cover to be removed.



6. Lift Cover vertically to remove from chassis.

