


Wasp Vacuum Cup Inspection

CAUTION

Heater block can be at HIGH TEMPERATURES DO NOT TOUCH by hand

Locate the Vacuum cups, positioned at the four corners of heater block, two at the front & two at rear.

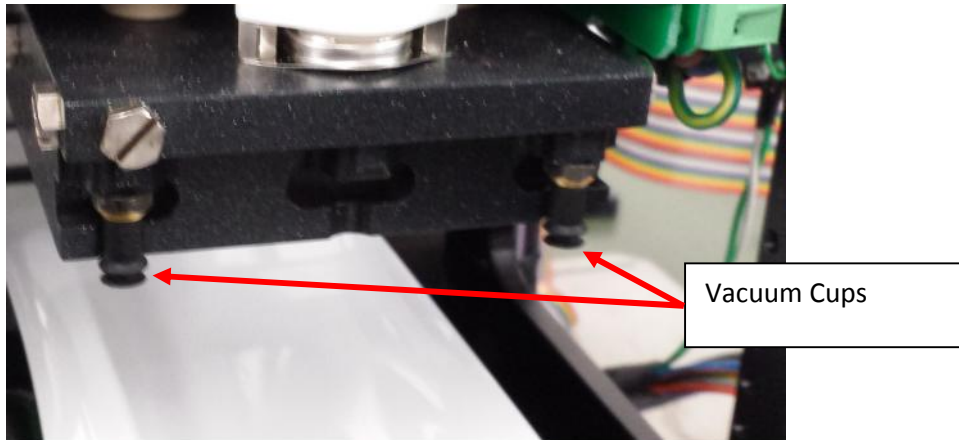

The vacuum cups must be of good condition to enable accurate and reliable foil/film pick up to continue with the sealing action of the Wasp plate sealer.

Each vacuum cup is a push fit onto the barbed fitting screwed into the heater block, ensure all fittings are complete and no damage has occurred.



Visually Inspect the Vacuum cup ensuring the 'sealing lip' of each vacuum cup is complete and not damaged.

