

Wasp Plate Sealer

Outer cover removal procedure

1. Disconnect power and Pneumatic services to Wasp Unit (Ensure Heater block is COOL)
2. Remove foil reel from rear of Unit.
3. Remove 3 screws and remove left hand side inspection cover.



4. Push Heater block down (**ensure it is cool before touching it**)



5. Remove the screen Bezel Assembly



6. To remove display, Tilt display backwards in opening and very carefully withdraw from cover.



Wasp

Outer cover removal procedure

7. Disconnect all connections from screen
 - a. Screen Ribbon Cable.
 - b. Two Pin Fan Lead
 - c. Thermocouple (two pin black connector Pulls off) Green wire is + positive



8. Remove display from cover.
9. Extreme care to be taken, do not separate the two Halves of touch screen!



10. Carefully lift cover from main chassis unit.



11. To replace cover, reverse this procedure.